



Procedure

1. Place the camera lens snug to the microscope eyepiece.
2. Orient the camera until you see the same image on your camera screen that you observed when looking through the microscope eyepiece.
3. Adjust the fine focus knob on the microscope as needed to ensure the image appears clear on your camera. See Figure 1.

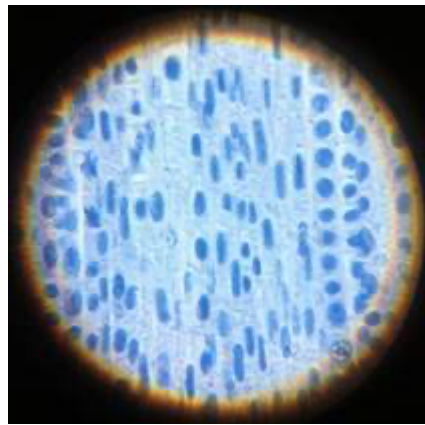


Figure 1. Onion root tip cells viewed on digital camera screen.

4. Take a photo of the focused slide.
5. Save the photo to your device in JPG, JPEG, PNG, or PDF format.
6. Label the image, if required, using PowerPoint, Google Slides, or a similar program.
7. Save the photo to your device in JPG, JPEG, PNG, or PDF format.
8. Upload the photo by selecting “Choose File” and selecting the image you saved. Then, click upload.